


Test System for Counting Bacteria and Total Coliform

DIRECTIONS FOR USE

1. Unscrew the cap and remove the paddle from the vial. Be careful not to touch either of the paddle's surfaces.
2. Immerse paddle into the sample for 2-3 seconds so that both surfaces are covered with liquid.
3. Allow excess fluid to drain from the paddle by touching the tip to an absorbant piece of paper (such as a paper towel).
4. Screw cap back and incubate SANI-CHECK BC in an upright position between 35 and 37° C ± for 24-48 hours.
5. Compare growth on paddle to the conversion chart below for a quantitative interpretation of the results.

Bacteria cfu / ml

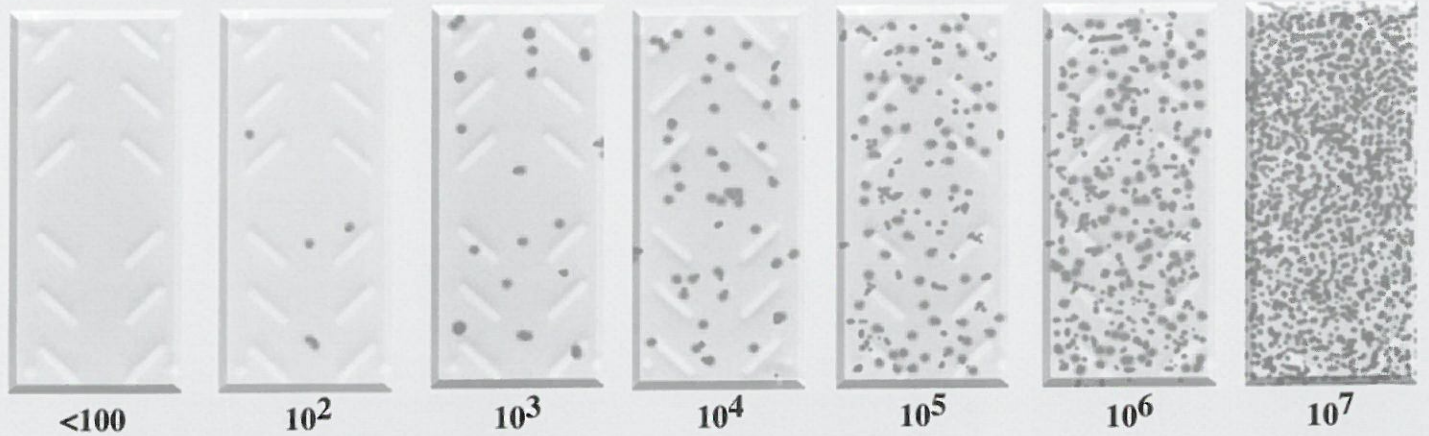

cfu/ml = colony forming units per milliliter

BIOSAN LABORATORIES, INC. offers technical service on matters relating to its products at no charge. Call toll free 800-253-6800 from within the U.S.A. and Canada, or 586-755-8970 from outside the U.S.A.



1950 Tobsal Court ■ Warren, MI 48091-1351 USA
Phone: 586.755.8970 ■ Toll Free: 800.253.6800 ■ Fax: 586.755.8978
www.biosan.com